



# Creating cDNA Libraries using Twist Library Preparation EF Kit 2.0 for ssRNA Virus Detection

For use with the Twist NGS Workflow

Most viral pathogens are single-stranded RNA including SARS-CoV-2, the virus behind the COVID-19 pandemic. This protocol describes how to convert RNA into cDNA libraries for downstream use in NGS Target Enrichment. The protocol begins with creating cDNA from extracted/purified RNA using commercially available reagents. The cDNA is then converted to Illumina TruSeq-compatible libraries using the Twist Library Preparation EF Kit 2.0 with Enzymatic Fragmentation and Unique Dual Indices (UDI). The resulting cDNA libraries are compatible with Twist Target Enrichment Protocols. Synthetic viral RNA controls, such as the Twist synthetic SARS-CoV-2 RNA controls, can be spiked into reference RNA samples to serve as a positive control.



## PROTOCOL COMPONENTS

Please read the product packaging and storage recommendations carefully for each kit and store components as recommended immediately upon arrival.

PART NUMBER	NAME	DESCRIPTION	STORAGE
See Appendix C	Twist Synthetic SARS-CoV-2 RNA Controls*	Synthetic SARS-CoV-2 RNA controls	-90 to -70°C
104206: 16 rxn 104207: 96 rxn	Twist Library Preparation EF Kit 2.0	Reagents for library construction	—
	Twist Library Preparation EF Kit 1, 2.0	Frag/AT Enzymes Frag/AT Buffer Ligation Master Mix Equinox Library Amp Mix (2x) P5/P7 Primers (10x)	-25 to -15°C
	Twist Library Preparation Kit 2	DNA Purification Beads	2 to 8°C
100401: 16 rxn 100573: 96 rxn	Twist Library Preparation Kit 2	DNA Purification Beads (as a Standalone Product, Bead Purification is also needed during cDNA Synthesis)	2 to 8°C
101307: 16 rxn 101308, 101309, 101310, 101311: 96 rxn	Twist Universal Adapter System - TruSeq Compatible	Twist Universal Adapters and Twist UDI Primers	-25 to -15°C

\*Optional as experimental control.

This product is for **research use only**.



## LEGAL

---

This document may contain references to other third-party resources such as sources of information, hardware or software, products, or services and/or web sites owned or licensed by third parties. Twist Bioscience does not control or take responsibility for any third-party resources, including, without limitation, the accuracy, reliability, copyright compliance, compatibility, performance, legality, or any other aspect of third-party resources. The inclusion of such resources in this document does not imply endorsement by Twist Bioscience of any third-party resources.

Certain processes described in this document may be subject to patent rights or licenses in local jurisdictions, including those owned or licensed by parties other than Twist Bioscience. Purchase of this product does not include a license to perform any such processes. Users of this product may, therefore, be required to obtain a patent license depending upon the particular application and country in which the product is used before performing such processes. Twist Bioscience, the various Twist logos, and the Twist NGS logo used herein are trademarks of Twist Bioscience Corporation.

All other trademarks are the property of their respective owners.

© 2021 Twist Bioscience Corporation. All rights reserved.

## INTENDED USE

---

This product is for research use only. This product is not intended for the diagnosis, prevention, or treatment of a disease or condition. Twist Bioscience assumes no liability regarding use of the product for applications in which it is not intended.



## TABLE OF CONTENTS

---

Materials Supplied by User	4
Equipment to be Supplied by User	5
General Notes and Precautions	6
Important Notes	6
Protocol Overview	7
Step 1: Dilution of Experimental RNA Samples and Synthetic Controls	8
• Dilute the Samples	8
Step 2: Execute cDNA Synthesis and Purify	9
• Prepare Primer Annealing Solution	9
• Prepare First Strand Synthesis Solution	10
• Prepare Second Strand Synthesis Solution	10
• Purify	10
Step 3: Perform DNA Fragmentation, End Repair, and dA-Tailing	12
• Prepare the Thermal Cycler, Samples, and Reagents	12
• Perform Fragmentation, End Repair, and dA-Tailing	13
Step 4: Ligate Twist Universal Adapters and Purify	14
• Ligate Twist Universal Adapters	14
• Purify	15
Step 5: PCR Amplify Using Twist UDI Primers, Purify, and Perform QC	16
• Prepare the Thermal Cycler and PCR Mix	16
• Perform PCR	17
• Purify	17
• Perform QC	18
Appendix A: Adapter Sequences and Pooling Guidelines	19
Appendix B: Twist Target Enrichment Protocols	21
Appendix C: Twist Synthetic SARS-CoV-2 Controls	23



# MATERIALS SUPPLIED BY USER

The following materials or their equivalent are required to generate libraries using the Twist Library Preparation EF Kit 2.0 for ssRNA kit.

PRODUCT	SUGGESTED SUPPLIER
ProtoScript II First Strand cDNA Synthesis Kit	New England Biolabs (E6560L, sufficient for 60 samples)
NEBNext Ultra II Non-Directional RNA Second Strand Synthesis Module	New England Biolabs (E6111)
Random Primer 6	New England Biolabs (S1230S)
Gene Expression Universal Reference RNA (Human)	Agilent Technologies (740000-41)
Ethanol (200 proof)	—
Molecular biology grade water	—
10 mM Tris-HCl pH 8 (optional)	—
Buffer EB (optional)	Qiagen
1.5-ml microcentrifuge tubes	VWR
Thin-walled PCR 0.2-ml strip-tubes	Eppendorf
96-well thermal cycling plates	VWR
96-well compatible magnetic plate	Alpaqua, Permagen Labware
Qubit RNA Broad Range Quantitation Assay	Thermo Fisher Scientific
Qubit dsDNA Broad Range Quantitation Assay	Thermo Fisher Scientific
Qubit dsDNA High Sensitivity Quantitation Assay	Thermo Fisher Scientific
Agilent DNA 7500 Kit	Agilent Technologies



# EQUIPMENT TO BE SUPPLIED BY USER

The following equipment or their equivalent are required to generate libraries using the Twist Library Preparation EF Kit 2.0 for ssRNA kit.

PRODUCT	SUGGESTED SUPPLIER
Pipettes and tips	—
Vortex mixer	—
Benchtop mini centrifuge for 0.2-ml tubes	—
Thermal cycler (96 well) with heated lid	—
Fluorometer (Qubit 3.0)	Thermo Fisher Scientific
2100 Bioanalyzer Agilent Technologies	Agilent Technologies



## GENERAL NOTES AND PRECAUTIONS

---

Wear appropriate protective equipment (lab coat, gloves, and protective glasses or goggles) at all times when performing this protocol.

For best results, read this document before performing the protocol and follow the instructions provided. Twist cannot guarantee performance if modifications are made to the protocol.

This library preparation method may yield more material than needed for target enrichment. Excess product can be stored at  $-20^{\circ}\text{C}$  for later use.

Test the compatibility of your thermal cycler and PCR tubes by incubating at  $95^{\circ}\text{C}$  for up to 5 minutes, to ensure the PCR tubes do not crack under heat and pressure. Adjust the tightness of the thermal cycler lid and/or use a spacer specific to the thermal cycler model.

## IMPORTANT NOTES

---

Store all RNA samples at  $-80^{\circ}\text{C}$  when not in use. Keep on ice at all times when in use.

Input RNA should be suspended in Molecular Biology Grade Water, 10 mM Tris-HCl pH 8.0, or Buffer EB. The recommended RNA input is 50 ng

Prior to library preparation, it is important to remove all cations and chelators as these may affect the initial fragmentation reaction.

For cDNA samples, correct input quantity is critical for achieving optimal yield and library fragment length.

To accommodate research samples with low viral titer, we recommend using 2.5X reactions of ProtoScript II First Strand cDNA Synthesis Kit compared to what NEB recommends using.

The recommended DNA input is 25 ng of purified cDNA.

Using higher cDNA input will increase library diversity and reduce duplication rate, but optimization of the following steps in library preparation may still be required to achieve optimal performance.

1. Incubation time for fragmentation (Step 3.1, page 12)
2. Amount of Twist Universal Adapter (Step 4.1, page 14)
3. Incubation time for ligation reaction (Step 4.4, page 14)

Measuring DNA concentration by absorbance at 260 nm is not recommended. Use the Thermo Fisher Scientific Qubit dsDNA Broad Range Quantitation Assay to accurately quantify input purified cDNA.

For technical support, contact [customersupport@twistbioscience.com](mailto:customersupport@twistbioscience.com).



## PROTOCOL OVERVIEW

This protocol begins with ssRNA and generates cDNA. This cDNA can then be amplified and indexed to create DNA libraries for subsequent target enrichment. The protocol features enzymatic fragmentation and Twist Universal Adapters with UDI primers. This protocol allows you to perform cDNA synthesis and library preparation (Steps 1–5) in 5 hours.

HUMAN GENOMIC RNA SAMPLES, HUMAN GENOMIC REFERENCE RNAS, AND TWIST SYNTHETIC CONTROLS		TIME
STEP 1	<b>Dilution of Experimental RNA Samples and Optional Synthetic Controls</b> Diluted Experimental RNA Samples and Synthetic Controls	10 minutes
STEP 2	<b>Execute cDNA Synthesis and Purify</b> cDNA material for Library Preparation	2.75 hours
cDNA, 50 NG STARTING DNA MATERIAL		
STEP 3	<b>Perform DNA Fragmentation, End Repair, and dA-Tailing</b> dA-Tailed DNA Fragments	1 hour
STEP 4	<b>Ligate Twist Universal Adapters and Purify</b> cDNA Libraries with Universal Adapters	1 hour
STEP 5	<b>PCR Amplify Using Twist UDI Primers, Purify, and Perform QC</b> Amplified Indexed Libraries	1 hour
STOPPING POINT		



## STEP 1

# DILUTION OF EXPERIMENTAL RNA SAMPLES AND SYNTHETIC CONTROLS

---

### Reagents Required

- Experimental RNA Samples: 50 ng RNA per sample
- Gene Expression Universal Reference RNA (Human)
- Twist Synthetic Controls
- Molecular Biology Grade Water

### Before You Begin

Thaw on ice:

- Experimental RNA Samples
- Gene Expression Universal Reference RNA (Human)
- Twist Synthetic Controls

## DILUTE THE SAMPLES

**1.1** Dilute the Experimental RNA Samples to 3.3 ng/μl in Molecular Biology Grade Water.

**1.2** Dilute the Gene Expression Universal Reference RNA (Human) to 3.85 ng/μl.

OPTIONAL: If using the Twist Synthetic RNA Control(s), we recommend diluting to the following concentrations:

- $0.5 \times 10^6$  copies/μl
- $0.5 \times 10^3$  copies/μl
- $0.5 \times 10^2$  copies/μl
- $0.5 \times 10^1$  copies/μl
- Negative Control (Molecular Biology Grade Water only)

**1.3** Add 2 μl of the diluted Twist Synthetic RNA Control to 13 μL of diluted Gene Expression Universal Reference RNA (Human) from step 1.2.

**PROCEED IMMEDIATELY TO STEP 2: EXECUTE cDNA SYNTHESIS AND PURIFY**





## STEP 2

## EXECUTE cDNA SYNTHESIS AND PURIFY

Generate cDNA from diluted Experimental RNA Samples and (optional) Synthetic Controls.

### Reagents Required

- Diluted Experimental RNA Samples and Synthetic Controls (from Step 1 - Samples are prepared in 1.1-1.3)
- Random Primer 6 (resuspended in water at 50  $\mu$ M)
- ProtoScript II First Strand cDNA Synthesis Kit
- NEBNext Ultra II Non-Directional RNA Second Strand Synthesis Module
- Ethanol
- Molecular Biology Grade Water
- From the Twist Library Preparation Kit 2:
  - DNA Purification Beads

### Before You Begin

- Thaw on ice:
  - ProtoScript II First Strand cDNA Synthesis Buffers and Reagents
  - NEBNext Ultra II Non-Directional RNA Second Strand Synthesis Buffers and Reagents
- Prepare 2 ml 80% ethanol for each sample (for use in Steps 2, 4 and 5 of the protocol).
- Equilibrate DNA Purification Beads to room temperature for at least 30 minutes (for use in Steps 2, 4 and 5 of the protocol).
- Resuspend the Random Primer 6 reagent with Molecular Biology Grade Water to 50  $\mu$ M.

## PREPARE PRIMER ANNEALING SOLUTION

### 2.1

Add the following volumes of reagents to each reaction tube (either a 0.2-ml thin walled PCR strip-tube or 96-well plate). Mix by flicking the tube(s).

REAGENT	VOLUME
Random Primer 6	5 $\mu$ l
Diluted Experimental RNA Samples or Synthetic Controls	15 $\mu$ l
Total	20 $\mu$ l

### 2.2

Pulse-spin the tube(s) and ensure there are minimal bubbles present.

### 2.3

Heat the Primer Annealing Solution to 95°C for 5 minutes in a thermal cycler with the lid at 105°C, then ramp down to 20°C. Immediately place on ice. Proceed to the next step.

## PREPARE FIRST STRAND SYNTHESIS SOLUTION

- 2.4** Add the following volumes of reagents to each tube of Primer Annealing Solution. Mix by flicking the tube(s).

REAGENT	VOLUME
ProtoScript II Reaction Mix	25 µl
ProtoScript II Enzyme Mix	5 µl

*Prepare a master mix for multiple reactions.*

- 2.5** Incubate the First Strand Synthesis Solution at the following conditions: 25°C for 5 minutes, 42°C for 1 hour, 80°C for 5 minutes, 4°C hold. Set the temperature of the heated lid to 105°C. When incubation is finished, proceed to the next step.

## PREPARE SECOND STRAND SYNTHESIS SOLUTION

- 2.6** Add the following volumes of reagents to each tube of First Strand Synthesis Solution. Mix by flicking the tube(s).

REAGENT	VOLUME
Molecular Biology Grade Water	18 µl
NEBNext Second Strand Synthesis Reaction Buffer	8 µl
NEBNext Second Strand Synthesis Enzyme Mix	4 µl

- 2.7** Incubate the Second Strand Synthesis Solution at 16°C for 1 hour in the thermal cycler, then move the samples to the bench top. Proceed to the Purify step.

 **IMPORTANT:** Turn off the heated lid or set to minimum temperature.

## PURIFY


- 2.8** Vortex the pre-equilibrated room temperature DNA Purification Beads until well mixed.
- 2.9** Add 96 µl of homogenized (1.2x) DNA Purification Beads to each Second Strand Synthesis Solution from Step 2.7. Mix well by vortexing.
- 2.10** Incubate the samples for 5 minutes at room temperature.
- 2.11** Place the samples on a magnetic plate for 1 minute or until the supernatant is clear.
- 2.12** The DNA Purification Beads form a pellet, leaving a clear supernatant. Without removing plate or tube(s) from the magnetic plate, remove and discard the supernatant.



- 2.13** Wash the bead pellet by gently adding 200  $\mu$ l freshly prepared 80% ethanol (do not disturb the pellet). Incubate for 1 minute, then remove and discard the ethanol.
- 2.14** Repeat the wash once, for a total of two washes, while keeping the sample(s) on the magnetic plate.
- 2.15** Carefully remove all remaining ethanol with a 10- $\mu$ l pipette, making sure not to disturb the bead pellet.  
**NOTE:** Before pipetting, the bead pellet may be briefly spun to collect ethanol at the bottom of the plate or tube and returned to the magnetic plate.
- 2.16** Air-dry the bead pellet on the magnetic plate for 5 minutes or until the bead pellet is dry. Do not overdry the bead pellet.
- 2.17** Remove the plate or tube(s) from the magnetic plate and add 27  $\mu$ l water to each sample. Mix by pipetting until homogenized.
- 2.18** Incubate at room temperature for 2 minutes.
- 2.19** Place the plate or tubes on a magnetic plate and let stand for 3 minutes or until the beads form a pellet.
- 2.20** Transfer 25  $\mu$ l of the clear supernatant containing the cDNA to a clean thin-walled PCR 0.2-ml strip-tube or well of a 96-well thermal cycling plate, making sure not to disturb the bead pellet.  
**OPTIONAL:** To perform a QC step, use the Thermo Fisher Scientific Qubit dsDNA High Sensitivity Quantification Assay.

**NOTE:** Final concentrations below 1 ng/ $\mu$ l may reflect inefficient sample preparation and can result in low library diversity after hybridization.

### PROCEED TO STEP 3: PERFORM DNA FRAGMENTATION, END REPAIR, AND dA-TAILING

 **STOPPING POINT:** If not immediately proceeding to Step 3, store the cDNA Pool at  $-20^{\circ}\text{C}$ .

## STEP 3 PERFORM DNA FRAGMENTATION, END REPAIR, AND dA-TAILING

Perform enzymatic fragmentation of input cDNA and subsequent end repair and dA-tailing to generate dA-tailed DNA fragments.

### Reagents Required

- cDNA (from Steps 2.20 above)
- Molecular Biology Grade Water
- Qubit dsDNA Broad Range Quantitation Assay (or equivalent)
- From the Twist Library Preparation EF Kit 1, 2.0:
  - Frag/AT Enzymes
  - Frag/AT Buffer

### Before You Begin

Thaw or place on ice:

- Molecular biology grade water
- cDNA
- Frag/AT Buffer
- Frag/AT Enzymes

## PREPARE THE THERMAL CYCLER, SAMPLES, AND REAGENTS

- 3.1** Program the thermal cycler with the following conditions. Use the Step 2 Incubation Time table to select conditions for fragmentation to achieve the desired insert size. Set the temperature of the heated lid to 105°C. Start the program to pre-chill the thermal cycler.

TEMPERATURE		TIME
STEP 1	4°C	HOLD
STEP 2	Use the table to the right to select the time and temperature for desired insert size	
STEP 3	65°C	30 minutes
STEP 4	4°C	HOLD

STEP 2 INCUBATION TIME*		
DESIRED INSERT SIZE (BP)	@30°C	@37°C
145–175	—	30 min
180–220	—	20 min
250–300	—	10 min
275–350	15 min	—
350–425	10 min	—

\*20 min at 37°C is recommended for Twist target enrichment applications. All recommendations are a starting point and should be optimized for each application.

- 3.2** Add 25 µl of each cDNA sample into either a thin-walled PCR 0.2-ml striptube or a well of a 96-well thermal cycling plate.

OPTIONAL: If QC step is performed after Step 2.20, dilute the cDNA samples to 1 ng/µl with water before aliquoting each cDNA sample out.

- 3.3** Pulse-spin to ensure all of the solution is at the bottom of the tube and place on ice.

## PERFORM FRAGMENTATION, END REPAIR, AND dA-TAILING

**3.4** Vortex the Frag/AT Buffer for 5 seconds. Pulse-spin to collect all liquid in the bottom of the tube.

**3.5** Invert Frag/AT Enzymes a minimum of 10 times to homogenize or briefly vortex to ensure complete mixing. Pulse-spin to collect all liquid in the bottom of the tube.

**3.6** Prepare an enzymatic fragmentation mix in a 1.5 ml microfuge tube on ice. Use the volumes listed below. Homogenize the mastermix with moderate vortexing for 5 seconds or pipetting a minimum of half the total volume up and down 10 times (avoid formation of bubbles).

REAGENT	VOLUME PER REACTION*
Molecular Biology Grade Water	15 µl
Frag/AT Buffer	4 µl
Frag/AT Enzymes	6 µl
Total	25 µl

*\*Prepare a master mix for multiple reactions.*

**3.7** Add 25 µl enzymatic fragmentation mix (from Step 3.6) to each 25 µl cDNA sample well or tube. Homogenize with moderate vortexing for 5 seconds or by pipetting a minimum of half the total volume up and down 10 times (avoid formation of bubbles). Cap the tube(s) or seal the plate and keep the reaction on ice.

**NOTE:** Complete mixing is critical to achieve consistent fragment lengths.

**3.8** Pulse-spin the sample plate or tube(s) and immediately transfer to the pre-chilled thermal cycler.

**3.9** Initiate steps 2 to 4 of the thermal cycler program (refer to the table in Step 3.1 above).

**NOTE:** While the thermal cycler program is running, prepare the reagents for Step 4: Ligate Twist Universal Adapters and Purify (see Before You Begin).

**3.10** When the thermal cycler program is complete and the sample block has returned to 4°C, remove the samples and place them on ice.

**PROCEED IMMEDIATELY TO STEP 4: LIGATE TWIST UNIVERSAL ADAPTERS AND PURIFY**

## STEP 4

# LIGATE TWIST UNIVERSAL ADAPTERS AND PURIFY

Ligate Twist Universal Adapters to the dA-tailed DNA fragments from Step 3 and purify to generate cDNA libraries ready for index introduction through amplification in Step 5.

### Reagents Required

- dA-tailed DNA fragments (from Step 3.10)
- 80% Ethanol (from Step 2)
- Equilibrated DNA Purification Beads (from Step 2)
- Molecular Biology Grade Water
- 10 mM Tris-HCl pH 8 or Buffer EB (optional, for elution)
- From the Twist Library Preparation EF Kit 1, 2.0:
  - Ligation Master Mix
- From the Twist Universal Adapter System:
  - Twist Universal Adapters
- From the Twist Library Preparation Kit 2:
  - DNA Purification Beads

### Before You Begin

- Thaw or place on ice:
  - Twist Universal Adapters (tube; utilized for all samples)
  - Ligation Master Mix
- Program a thermal cycler to incubate samples at 20°C with the heated lid set to minimum temperature or turned off. Start the program so that the cycler has reached 20°C when the samples are done being prepared.

## LIGATE TWIST UNIVERSAL ADAPTERS

**4.1** Add 2.5 µl of Twist Universal Adapters into each sample well or tube containing the dA-tailed DNA fragments from Step 3.10. Mix gently by pipetting and keep on ice.

**4.2** Invert the Ligation Master Mix a minimum of 10 times until homogenized and place on ice.

**NOTE:** Do not vortex the Ligation Master Mix.

**4.3** Add 2.5 µl of Molecular Biology Grade Water and 20 µl of Ligation Master Mix to each sample from Step 4.1. Pipette a minimum of half the total volume up and down 10 times to ensure complete mixing. Seal or cap the sample plate or tube(s) and pulse-spin to ensure all solution is at the bottom of the tube.

**4.4** Incubate the ligation reaction at 20°C for 15 minutes in the thermal cycler, then move the samples to the bench top. Proceed to the Purify step.

**⚠ IMPORTANT:** Turn off the heated lid or set to minimum temperature.

**NOTE:** While the thermal cycler program is running, prepare the reagents for Step 5: PCR Amplify Using Twist UDI Primers, Purify, and Perform QC (see Before You Begin).

**PURIFY**

- 4.5** Vortex the pre-equilibrated room temperature DNA Purification Beads until well mixed.
- 4.6** Add 60  $\mu$ l of homogenized (0.8x) DNA Purification Beads to each ligation sample from Step 4.4. Mix well by vortexing.
- 4.7** Incubate the samples for 5 minutes at room temperature
- 4.8** Place the samples on a magnetic plate for 1 minute or until the supernatant is clear.
- 4.9** The DNA Purification Beads form a pellet, leaving a clear supernatant. Without removing plate or tube(s) from the magnetic plate, remove and discard the supernatant.
- 4.10** Wash the bead pellet by gently adding 200  $\mu$ l freshly prepared 80% ethanol (do not disturb the pellet). Incubate for 1 minute, then remove and discard the ethanol.
- 4.11** Repeat the wash once, for a total of two washes, while keeping the sample(s) on the magnetic plate.
- 4.12** Carefully remove all remaining ethanol with a 10- $\mu$ l pipette, making sure not to disturb the bead pellet.  
**NOTE:** Before pipetting, the bead pellet may be briefly spun to collect ethanol at the bottom of the plate or tube and returned to the magnetic plate.
- 4.13** Air-dry the bead pellet on the magnetic plate for 5 minutes or until the bead pellet is dry. Do not overdry the bead pellet.
- 4.14** Remove the plate or tube(s) from the magnetic plate and add 17  $\mu$ l water to each sample. Mix by pipetting until homogenized.  
**NOTE:** 10 mM Tris-HCl pH 8 or Buffer EB may also be utilized for elution.
- 4.15** Incubate at room temperature for 2 minutes.
- 4.16** Place the plate or tubes on a magnetic plate and let stand for 3 minutes or until the beads form a pellet.
- 4.17** Transfer 15  $\mu$ l of the clear supernatant containing the ligated and indexed libraries to a clean thin-walled PCR 0.2 ml strip-tube or well of a 96-well thermal cycling plate, making sure not to disturb the bead pellet.

**PROCEED TO STEP 5: PCR AMPLIFY USING TWIST UDI PRIMERS, PURIFY, AND PERFORM QC**

## STEP 5

# PCR AMPLIFY USING TWIST UDI PRIMERS, PURIFY, AND PERFORM QC

Complete the protocol by amplifying the adapted cDNA libraries with Twist UDI Primers, purifying them, and performing QC analysis.

### Reagents Required

- Ligated, adapted libraries (from Step 4.17)
- 80% Ethanol (from Step 4)
- Equilibrated DNA Purification Beads (from Step 4)
- Molecular biology grade water
- 10 mM Tris-HCl pH 8 or Buffer EB (optional, for elution)
- From the Twist Library Preparation EF Kit 1, 2.0:
  - Equinox Library Amp Mix (2x)
- From the Twist Universal Adapter System:
  - Twist UDI Primers

**⚠ IMPORTANT:** Use of P5/P7 Primers (10x) tubes 104127, 103842 contained in the Twist Library Preparation EF Kit 1, 2.0 is not required. Using these primers with the Twist Universal Adapter System will result in a failed PCR amplification.

### Before You Begin

Thaw or place on ice:

- Equinox Library Amp Mix (2x)
- Twist UDI Primers (plate with single-use primers)

## PREPARE THE THERMAL CYCLER AND PCR MIX

### 5.1

Program a thermal cycler with the following conditions. Set the temperature of the heated lid to 105°C.

STEP	TEMPERATURE	TIME	NUMBER CYCLES
1 Initialization	98°C	45 seconds	1
2 Denaturation  Annealing  Extension	98°C	15 seconds	8-12*
	60°C	30 seconds	
	72°C	30 seconds	
3 Final Extension	72°C	1 minute	1
4 Final Hold	4°C	HOLD	—

\*Number of cycles will depend on the integrity and purity of starting cDNA.



**PERFORM PCR**

**5.2** Add 10 µl of Twist UDI Primer from the provided 96-well plate to each of the cDNA libraries from Step 4.17 and mix well by gentle pipetting.

**NOTE:** For index selection and pooling guidelines for downstream target enrichment and sequencing, refer to the Appendix.

**5.3** Add 25 µl of Equinox Library Amp Mix (2x) to the cDNA libraries from Step 5.2 and mix well by gentle pipetting.

**NOTE:** Invert Equinox Library Amp Mix (2x) 5 times before use. Do not vortex.

**5.4** Pulse-spin sample plate or tube and immediately transfer to the thermal cycler. Start the program.

**5.5** Remove the sample(s) from the block when the thermal cycler program is complete. Proceed to purification.

**PURIFY**

**5.6** Vortex the pre-equilibrated DNA Purification Beads until mixed.

**5.7** Add 50 µl (1x) of homogenized DNA Purification Beads to each ligation sample from Step 5.5. Mix well by vortexing.

**5.8** Incubate the samples for 5 minutes at room temperature.

**5.9** Place the samples on a magnetic plate for 1 minute.

**5.10** The DNA Purification Beads form a pellet, leaving a clear supernatant. With plate or tubes still on the magnetic plate, remove and discard the supernatant.

**5.11** Gently wash the bead pellet by adding 200 µl of freshly prepared 80% ethanol (do not disturb the pellet), incubate for 1 minute, then remove and discard the ethanol.

**5.12** Repeat this wash once, for a total of two washes, while keeping the samples on the magnetic plate.

**5.13** Carefully remove all remaining ethanol with a 10 µl pipet, making sure not to disturb the pellet.

**5.14** Air-dry the pellet on the magnetic plate for 5–10 minutes or until it is dry. Do not overdry.

**5.15** Remove the plate or tubes from the magnetic plate and add 22 µl of water, 10 mM Tris-HCl pH 8, or Buffer EB to each sample. Mix by pipetting until homogenized.


**5.16** Incubate at room temperature for 2 minutes.

**5.17** Place the plate or tubes on a magnetic plate and let stand for 3 minutes or until the beads form a pellet.

**5.18** Transfer 20 µl of the clear supernatant containing the Amplified Indexed Libraries to a clean thin-walled PCR 0.2 ml strip-tube or well of a 96-well thermal cycling plate, making sure not to disturb the pellet.

## PERFORM QC

**5.19** Quantify and validate the size range of each library using the Thermo Fisher Scientific Qubit dsDNA Broad Range Quantitation Assay and Agilent DNA 7500 Assay. Average fragment length should be 250–450 bp using a range setting of 150–1000 bp.

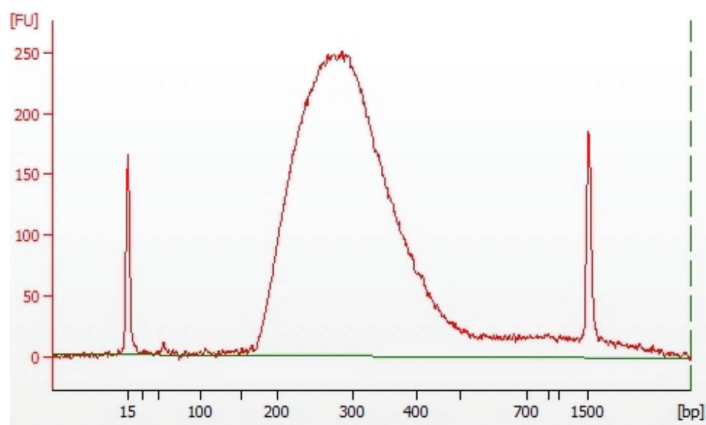

*Electropherogram generated by an Agilent 7500 DNA analysis of cDNA library samples that were prepared as described. Note the single prominent peak.*

**NOTES:** If the average fragment length is not within the desired range, optimize the fragmentation conditions used in Step 3.1. Test a range of fragmentation times in 3-5 minute increments around the target size from the Step 2 Incubation Time table. Increase time to produce shorter fragments and decrease time to produce longer fragments.

**STOPPING POINT:** If not immediately proceeding to a Twist Target Enrichment System, store the amplified indexed libraries at  $-20^{\circ}\text{C}$ .

## END OF WORKFLOW



# APPENDIX A: ADAPTER SEQUENCES AND POOLING GUIDELINES

## UDI SEQUENCES

For a complete guide of the Twist UDI sequences please refer to PDF document DOC-001129 or Excel file DOC-001130. Both files are available for download at [twistbioscience.com/resources](https://www.twistbioscience.com/resources).

## POOLING GUIDELINES

Twist UDI primers are base balanced by column for next generation sequencing. When pooling unique dual-indexed libraries for 8-plex hybridization, it is recommended that libraries be selected from a single column. Multiple columns may be selected in any desired combination across a single plate or multiple plates for sequencing.

**Table 1. Twist UDI primer plate layouts and pooling guidelines.**  
Twist Universal Adapter System: TruSeq Compatible, 16 Samples (101307)

	1	2	3	4	5	6	7	8	9	10	11	12
A	1	9										
B	2	10										
C	3	11										
D	4	12										
E	5	13										
F	6	14										
G	7	15										
H	8	16										

\*PLEASE NOTE: The indexes in the 16-sample plate are not the same for 96 samples (Plate A).



## APPENDIX A: ADAPTER SEQUENCES AND POOLING GUIDELINES

Twist Universal Adapter System: TruSeq Compatible, 96 Samples, Plates A to D (101308, 101309, 101310, 101311)

**Plate A.**

	1	2	3	4	5	6	7	8	9	10	11	12
A	1	9	17	25	33	41	49	57	65	73	81	89
B	2	10	18	26	34	42	50	58	66	74	82	90
C	3	11	19	27	35	43	51	59	67	75	83	91
D	4	12	20	28	36	44	52	60	68	76	84	92
E	5	13	21	29	37	45	53	61	69	77	85	93
F	6	14	22	30	38	46	54	62	70	78	86	94
G	7	15	23	31	39	47	55	63	71	79	87	95
H	8	16	24	32	40	48	56	64	72	80	88	96

**Plate B.**

	1	2	3	4	5	6	7	8	9	10	11	12
A	97	105	113	121	129	137	145	153	161	169	177	185
B	98	106	114	122	130	138	146	154	162	170	178	186
C	99	107	115	123	131	139	147	155	163	171	179	187
D	100	108	116	124	132	140	148	156	164	172	180	188
E	101	109	117	125	133	141	149	157	165	173	181	189
F	102	110	118	126	134	142	150	158	166	174	182	190
G	103	111	119	127	135	143	151	159	167	175	183	191
H	104	112	120	128	136	144	152	160	168	176	184	192

**Plate C.**

	1	2	3	4	5	6	7	8	9	10	11	12
A	193	201	209	217	225	233	241	249	257	265	273	281
B	194	202	210	218	226	234	242	250	258	266	274	282
C	195	203	211	219	227	235	243	251	259	267	275	283
D	196	204	212	220	228	236	244	252	260	268	276	284
E	197	205	213	221	229	237	245	253	261	269	277	285
F	198	206	214	222	230	238	246	254	262	270	278	286
G	199	207	215	223	231	239	247	255	263	271	279	287
H	200	208	216	224	232	240	248	256	264	272	280	288

**Plate D.**

	1	2	3	4	5	6	7	8	9	10	11	12
A	289	297	305	313	321	329	337	345	353	361	369	377
B	290	298	306	314	322	330	338	346	354	362	370	378
C	291	299	307	315	323	331	339	347	355	363	371	379
D	292	300	308	316	324	332	340	348	356	364	372	380
E	293	301	309	317	325	333	341	349	357	365	373	381
F	294	302	310	318	326	334	342	350	358	366	374	382
G	295	303	311	319	327	335	343	351	359	367	375	383
H	296	304	312	320	328	336	344	352	360	368	376	384

**END OF APPENDIX A**



## APPENDIX B: TWIST TARGET ENRICHMENT PROTOCOLS

The Twist Standard Target Enrichment Protocol is compatible with this workflow. It generates enriched TruSeq DNA libraries for sequencing on Illumina next-generation sequencing (NGS) systems using a 16-hour hybridization time in a two-day target enrichment workflow. This protocol can be found at [twistbioscience.com/resources](https://twistbioscience.com/resources). When using the Twist SARS-CoV-2 Research Panel in conjunction with this target enrichment workflow, there are no modifications.

PART NUMBER	NAME	DESCRIPTION	STORAGE
102016: 2 rxn 102017: 12 rxn 102018: 96 rxn	Twist SARS-CoV-2 Research Panel	Custom DNA Panels for SARS-CoV-2 Viral Detection	-25 to -15°C
100856: 2 rxn 100578: 12 rxn 100767: 96 rxn	Twist Universal Blockers	For the prevention of nonspecific capture: Universal Blockers Blocker Solution	-25 to -15°C
101262: 2 rxn 100983: 12 rxn 100984: 96 rxn	Twist Binding and Purification Beads	For target enrichment and purification: Streptavidin Binding Beads DNA Purification Beads	2 to 8°C
101279: 2 rxn 101025: 12 rxn 101026: 96 rxn	Twist Hybridization and Wash Kit (2 Boxes)	For target enrichment with standard hybridization:	-25 to -15°C  2 to 8°C
	Twist Hybridization Reagents (Box 1 of 2)	Hybridization Mix Hybridization Enhancer Amplification Primers	
	Twist Wash Buffers (Box 2 of 2)	Binding Buffer Wash Buffer 1 Wash Buffer 2	



## APPENDIX B: TWIST TARGET ENRICHMENT PROTOCOLS

The Twist Fast Hybridization Target Enrichment system is also compatible for viral detection use. This workflow uses hybridization times between 30 minutes and 2 hours, and is a one-day target enrichment workflow. This protocol can be found at [twistbioscience.com/resources](https://twistbioscience.com/resources). When using the Twist SARS-CoV-2 Research Panel in conjunction with this target enrichment workflow, the Fast Wash Buffer 1 temperature should be set at 68°C.

PART NUMBER	NAME	DESCRIPTION	STORAGE
102016: 2 rxn 102017: 12 rxn 102018: 96 rxn	Twist SARS-CoV-2 Research Panel	Custom DNA Panels for SARS-CoV-2 Viral Detection	-25 to -15°C
100856: 2 rxn 100578: 12 rxn 100767: 96 rxn	Twist Universal Blockers	For the prevention of nonspecific capture: Universal Blockers Blocker Solution	-25 to -15°C
101262: 2 rxn 100983: 12 rxn 100984: 96 rxn	Twist Binding and Purification Beads	For target enrichment and purification: Streptavidin Binding Beads DNA Purification Beads	2 to 8°C
101278: 2 rxn 101174: 12 rxn 101175: 96 rxn	Twist Fast Hybridization and Wash Kit (2 Boxes)	For target enrichment with fast hybridization:	-25 to -15°C  2 to 8°C
	Twist Fast Hybridization Reagents (Box 1 of 2)	Fast Hybridization Mix Hybridization Enhancer Amplification Primers	
	Twist Fast Wash Buffers (Box 2 of 2)	Fast Binding Buffer Fast Wash Buffer 1 Wash Buffer 2	

END OF APPENDIX B



## APPENDIX C: TWIST SYNTHETIC SARS-COV-2 CONTROLS

---

Each viral control is made synthetically with Twist's silicon platform and *in vitro* transcription. Each control is broken up between six unique 5 kb fragments, spanning the viral genome. For a full list of available viral controls, please visit: [twistbioscience.com/resources/product-sheet/Twist-Synthetic-Infectious-Disease-Controls-Product-List](https://twistbioscience.com/resources/product-sheet/Twist-Synthetic-Infectious-Disease-Controls-Product-List)

### END OF APPENDIX C

LAST REVISED: November 1, 2022

REVISION	DATE	DESCRIPTION
6.0	Nov 1, 2022	<ul style="list-style-type: none"><li>• Language changes to clarify reagent handling, required inputs, and steps</li><li>• Updated insert size optimization guidelines</li><li>• Removed table in Appendix C and provided website link with updated synthetic controls list</li></ul>
5.0	Jan 14, 2022	<ul style="list-style-type: none"><li>• Minor language update to clarify workflow steps.</li></ul>
4.0	Oct 28, 2021	<ul style="list-style-type: none"><li>• Protocol updated from SARS-CoV-2 workflow to a more general total nucleic acid protocol</li><li>• Minor language change to recommend optimizing fragmentation conditions depending on sample type, application and concentration</li></ul>
3.0	Sep 24, 2021	<ul style="list-style-type: none"><li>• Protocol updated with Twist Library Preparation EF Kit 2.0 workflow</li><li>• Minor updates to clarify workflow steps</li></ul>