



LIBRARY PREPARATION PROTOCOL

Enzymatic Fragmentation and Twist UMI Adapter System

For use with the Twist NGS Workflow

This Twist Library Preparation Kit provides the reagents needed to prepare genomic DNA (gDNA) libraries using enzymatic gDNA fragmentation and unique molecular identifiers (UMIs). The Twist UMI Adapter system consists of Twist UMI Adapters and Twist Unique Dual Indexed (UDI) Primers. This manual details the steps for generating the amplified, indexed libraries with UMI tags for downstream target enrichment and sequencing on Illumina next-generation sequencing (NGS) systems. This library preparation protocol is optimized for use with Twist Target Enrichment Kits and should only be performed with reagents specified or their equivalents.



Twist NGS workflow. The complete NGS workflow takes you from sample preparation to NGS sequencing and data analysis. A component of this workflow, the Twist Library Preparation Protocol works in conjunction with the other component protocols.

PROTOCOL COMPONENTS

Please read the product packaging and storage recommendations carefully for each kit and store components as recommended immediately upon arrival.

CATALOG #	NAME	DESCRIPTION	STORAGE
104174: 16 rxn 104175: 96 rxn	Twist Library Preparation EF Kit with Amp Mix	Reagents for library construction	—
	Twist Library Preparation EF Kit 1	· 5x Fragmentation Enzyme · 10x Fragmentation Buffer · DNA Ligation Mix · DNA Ligation Buffer · Amplification Primers, ILMN (Tubes 100220, 100583 are not required when used with UMI adapters)	–20°C
	Twist Library Preparation Kit 2	DNA Purification Beads	2–8°C
	Equinox Library Amp Mix	Equinox Library Amp Mix (2x)	–20°C
105040: 16 rxn 105041, 105042, 105043, 105044: 96 rxn	Twist UMI Adapter System - TruSeq Compatible	Twist UMI Adapters and Twist UDI Primers provide UMI labeling and unique dual-indexed combinations with 1 reaction per index pair	–20°C

This product is for **research use only**.



LEGAL

This document may contain references to other third-party resources such as sources of information, hardware or software, products, or services and/or web sites owned or licensed by third parties. Twist Bioscience does not control or take responsibility for any third-party resources, including, without limitation, the accuracy, reliability, copyright compliance, compatibility, performance, legality, or any other aspect of third-party resources. The inclusion of such resources in this document does not imply endorsement by Twist Bioscience of any third-party resources.

Certain processes described in this document may be subject to patent rights or licenses in local jurisdictions, including those owned or licensed by parties other than Twist Bioscience. Purchase of this product does not include a license to perform any such processes. Users of this product may, therefore, be required to obtain a patent license depending upon the particular application and country in which the product is used before performing such processes.

Twist Bioscience, the various Twist logos, and the Twist NGS logo used herein are trademarks of Twist Bioscience Corporation. All other trademarks are the property of their respective owners.

©2023 Twist Bioscience Corporation. All rights reserved.

INTENDED USE

This product is for research use only. This product is not intended for the diagnosis, prevention, or treatment of a disease or condition. Twist Bioscience assumes no liability regarding use of the product for applications in which it is not intended.



TABLE OF CONTENTS

Twist Library Preparation Protocol	1
Materials Supplied by User	4
General Notes and Precautions	5
Guidelines for gDNA Samples	5
Protocol Overview	6
Step 1: Perform DNA Fragmentation, End Repair, and dA-Tailing	7
• Prepare the Thermal Cycler, Samples, and Reagents	7
• Perform Fragmentation, End Repair, and dA-Tailing	8
Step 2: Ligate Twist UMI Adapters and Purify	9
• Ligate Twist UMI Adapters	9
• Purify	10
Step 3: PCR Amplify Using Twist UDI Primers, Purify, and Perform QC	11
• Prepare the Thermal Cycler and PCR Mix	11
• Perform PCR	12
• Purify	12
• Perform QC	13
Appendix: UDI Sequences and Pooling Guidelines	14



MATERIALS SUPPLIED BY USER

The following materials or their equivalent are required to generate libraries using the Twist Library Preparation Kit with Enzymatic Fragmentation and Twist UMI Adapter System.

PRODUCT	SUGGESTED SUPPLIER
REAGENTS AND CONSUMABLES	
Ethanol (200 proof)	—
Molecular biology grade water	—
10 mM Tris-HCl pH 8	—
Buffer EB	Qiagen
1.5-ml microcentrifuge tubes	VWR
Thin-walled PCR 0.2-ml strip-tubes	Eppendorf
96-well thermal cycling plates	VWR
1.5-ml compatible magnetic stand	Beckman Coulter
96-well compatible magnetic plate	Alpaqua, Permagen Labware
Qubit dsDNA Broad Range Quantitation Assay	Thermo Fisher Scientific
Agilent DNA 7500 Kit	Agilent Technologies
EQUIPMENT	
Pipettes and tips	—
Vortex mixer	—
Benchtop mini centrifuge for 0.2-ml tubes	—
Thermomixer for 1.5-ml tubes	Eppendorf
Thermal cycler (96 well) with heated lid	—
Fluorometer (Qubit 3.0)	Thermo Fisher Scientific
2100 Bioanalyzer	Agilent Technologies



GENERAL NOTES AND PRECAUTIONS

Wear appropriate protective equipment (lab coat, gloves, and protective glasses or goggles) at all times when performing this protocol.

For best results, read this document before performing the protocol, and follow the instructions provided. Twist cannot guarantee the performance of the Twist Library Preparation Kit using Enzymatic Fragmentation and the Twist UMI Adapter System if modifications are made to the protocol.

This library preparation method may yield more material than needed for target enrichment. Excess product can be stored at -20°C for later use.

Do NOT mix or combine the same reagents from the different lots.

Test the compatibility of your thermal cycler and PCR tubes by incubating at 95°C for up to 5 minutes to ensure the PCR tubes do not crack under heat and pressure. Adjust the tightness of the thermal cycler lid and/or use a spacer specific to the thermal cycler model.

GUIDELINES FOR gDNA SAMPLES

- Input DNA should be suspended in Molecular Biology Grade Water, 10 mM Tris-HCl pH 8.0, or Buffer EB.
- It is important to remove all cations and chelators from the starting gDNA sample. The presence of cations and chelators may affect the initial fragmentation reaction.
- For genomic (gDNA) samples, correct input quantity is critical for achieving optimal yield and library fragment length.
- The recommended DNA input is 50 ng of purified gDNA.
- Using higher or lower mass input may require optimization of the following steps in library preparation to achieve optimal performance.
 - Incubation time for fragmentation at 32°C (Step 1.1, page 7)
 - Amount of Twist UMI Adapter (Step 2.1, page 9)
 - Incubation time for ligation reaction (Step 2.5, page 10)
- Use the Thermo Fisher Scientific Qubit dsDNA Broad Range Quantitation Assay to accurately quantify input purified gDNA.
- Measuring DNA concentration by absorbance at 260 nm is not recommended.
- For technical support, contact customersupport@twistbioscience.com.



PROTOCOL OVERVIEW

This protocol begins with genomic DNA (gDNA) and generates amplified, indexed libraries for subsequent target enrichment. It features enzymatic fragmentation and Twist UMI Adapters with UDI primers. This protocol allows you to perform gDNA library preparation (Steps 1–3) in 3 hours.

ENZYMATIC FRAGMENTATION WITH UMI ADAPTERS AND UDI PRIMERS (GENOMIC DNA, 50 NG STARTING DNA MATERIAL)		TIME
STEP 1	Perform DNA Fragmentation, End Repair, and dA-tailing dA-tailed DNA fragments	1 hour
STEP 2	Ligate Twist UMI Adapters and Purify gDNA libraries ready for indexing	1 hour
STEP 3	PCR Amplify Using Twist UDI Primers, Purify, and Perform QC Amplified indexed libraries	1 hour



STEP 1 PERFORM DNA FRAGMENTATION, END REPAIR, AND dA-TAILING

Perform enzymatic fragmentation of input gDNA and subsequent end repair and dA-tailing to generate dA-tailed DNA fragments.

Reagents Required

- Genomic DNA (gDNA): 50 ng per sample
- Molecular biology grade water (chilled)
- Optional 10 mM Tris-HCl pH 8 or Buffer EB
- Qubit dsDNA Broad Range Quantitation Assay (or equivalent)
- From the Twist Library Preparation EF Kit 1:
 - 5x Fragmentation Enzyme
 - 10x Fragmentation Buffer

Before You Begin

- Thaw 5x Fragmentation Enzyme and gDNA samples on ice, then mix by flicking the tube with a finger.
- Thaw 10x Fragmentation Buffer on ice, then mix by pulse vortexing for 2 seconds.

PREPARE THE THERMAL CYCLER, SAMPLES, AND REAGENTS

- 1.1** Program the thermal cycler with the following conditions. Set the temperature of the heated lid to 70°C. Start the program to pre-chill the thermal cycler.

	TEMPERATURE	TIME
STEP 1	4°C	HOLD
STEP 2	32°C	22 minutes (adjustment may be required based upon DNA mass input and/or quality)
STEP 3	65°C	30 minutes
STEP 4	4°C	HOLD

- 1.2** Use the Qubit dsDNA Broad Range Quantitation Assay to determine the concentration of your gDNA samples.

- 1.3** Dilute the gDNA samples to 5 ng/μl with water, 10 mM Tris-HCl pH 8, or Buffer EB. Mix well with gentle pipetting.

- 1.4** Add 10 μl of each diluted gDNA sample (50 ng total gDNA) into a thin-walled PCR 0.2-ml strip-tube or well of a 96-well thermal cycling plate, and place on ice.

- 1.5** Pulse-spin to ensure all of the solution is at the bottom of the tube.

**PERFORM FRAGMENTATION, END REPAIR, AND dA-TAILING**

- 1.6** Prepare an enzymatic fragmentation master mix in a 1.5-ml microcentrifuge tube on ice. Use the volumes listed below. Mix thoroughly by gentle pipetting.

REAGENT	VOLUME PER REACTION*
Water (chilled)	25 µl
10x Fragmentation Buffer	5 µl
5x Fragmentation Enzyme	10 µl
Total	40 µl

**Prepare a master mix for multiple reactions.*

- 1.7** Add 40 µl enzymatic fragmentation master mix (from Step 1.6) to each 10 µl gDNA sample well or tube and mix well by gentle pipetting. Cap the tube and keep the reaction on ice.

- 1.8** Pulse-spin the sample plate or tubes and immediately transfer to the pre-chilled thermal cycler.

- 1.9** Proceed to steps 2–4 of the thermal cycler program (32°C step of the thermocycler program in Step 1.1 above).

NOTE: While the thermal cycler program is running, prepare the reagents for Step 2: Ligate Twist UMI Adapters and Purify (see Before You Begin).

- 1.10** When the thermal cycler program is complete and the sample block has returned to 4°C, remove the samples from the block and place on ice.

PROCEED IMMEDIATELY TO STEP 2: LIGATE TWIST UMI ADAPTERS AND PURIFY

STEP 2

LIGATE TWIST UMI ADAPTERS AND PURIFY

Ligate Twist UMI Adapters to the dA-tailed DNA fragments from Step 1 and purify to generate gDNA libraries ready for index introduction through amplification in Step 3.

Reagents Required

- dA-tailed DNA fragments (from Step 1.10)
- Ethanol
- Molecular biology grade water (chilled)
- Optional 10 mM Tris-HCl pH 8 or Buffer EB
- From the Twist Library Preparation EF Kit 1:
 - DNA Ligation Mix
 - DNA Ligation Buffer
- From the Twist UMI Adapter System:
 - Twist UMI Adapters
- From the Twist Library Preparation Kit 2:
 - DNA Purification Beads

Before You Begin

- Thaw on ice:
 - Twist UMI Adapters (tube; utilized for all samples)
 - DNA Ligation Mix
 - DNA Ligation Buffer
- Prepare 1 ml 80% ethanol for each sample (for use in Steps 2 and 3 of the protocol).
- Equilibrate DNA Purification Beads to room temperature for at least 30 minutes (for use in Steps 2 and 3 of the protocol).
- Program a thermal cycler to incubate the samples at 20°C with the heated lid set to minimum temperature or turned off. Start the program so that the cycler is at 20°C when the samples are prepared.

LIGATE TWIST UMI ADAPTERS

- 2.1** Add 3 µl Twist UMI Adapters into each sample well or tube containing the dA-tailed DNA fragments for Step 1. Mix gently by pipetting and keep on ice.
- 2.2** Prepare the ligation master mix in a 1.5-ml microcentrifuge tube on ice as indicated below. Mix well by gentle pipetting.

REAGENT	VOLUME PER REACTION*
Water (chilled)	17 µl
DNA Ligation Buffer	20 µl
DNA Ligation Mix	10 µl
Total	47 µl

*Prepare a master mix for multiple reactions.



- 2.3** Add 47 µl of the ligation master mix to the sample from Step 2.1 and mix well by gentle pipetting.
- 2.4** Seal or cap the tubes and pulse-spin to ensure all solution is at the bottom of the tube.
- 2.5** Incubate the ligation reaction at 20°C for 15 minutes in the thermal cycler, then move the samples to the bench top. Proceed to the Purify step.

⚠ IMPORTANT: Turn off the heated lid or set to minimum temperature.

NOTE: While the thermal cycler program is running, prepare the reagents for Step 3 (see Step 3: PCR Amplify Using Twist UDI Primers, Purify, and Perform QC).

PURIFY

- 2.6** Vortex the pre-equilibrated DNA Purification Beads until well mixed.
- 2.7** Add 80 µl (0.8x) of homogenized DNA Purification Beads to each ligation sample from Step 2.5. Mix well by vortexing.
- 2.8** Incubate the samples for 5 minutes at room temperature.
- 2.9** Place the samples on a magnetic plate for 1 minute or until the supernatant is clear.
- 2.10** The DNA Purification Beads form a pellet, leaving a clear supernatant. Without removing plate or tubes from the magnetic plate, remove and discard the supernatant.
- 2.11** Wash the bead pellet by gently adding 200 µl freshly prepared 80% ethanol (do not disturb the pellet). Incubate for 1 minute, then remove and discard the ethanol.
- 2.12** Repeat the wash once, for a total of two washes, while keeping the samples on the magnetic plate.
- 2.13** Carefully remove all remaining ethanol with a 10-µl pipet, making sure not to disturb the bead pellet.
- NOTE:** Before pipetting, the bead pellet may be briefly spun to collect ethanol at the bottom of the plate or tube and returned to the magnetic plate.
- 2.14** Air-dry the bead pellet on the magnetic plate for 5 minutes or until the bead pellet is dry. Do not overdry the bead pellet.
- 2.15** Remove the plate or tubes from the magnetic plate and add 17 µl water, 10 mM Tris-HCl pH 8, or Buffer EB to each sample. Mix by pipetting until homogenized.
- 2.16** Incubate at room temperature for 2 minutes.
- 2.17** Place the plate or tubes on a magnetic plate and let stand for 3 minutes or until the beads form a pellet.
- 2.18** Transfer 15 µl of the clear supernatant containing the ligated and indexed libraries to a clean thin-walled PCR 0.2-ml strip-tube or well of a 96-well thermal cycling plate, making sure not to disturb the bead pellet.

STEP 3

PCR AMPLIFY USING TWIST UDI PRIMERS, PURIFY, AND PERFORM QC

Amplify the adapted gDNA libraries with Twist UDI Primers, purify them, and perform quality control (QC) analysis to complete the protocol.

Reagents Required

- Ligated, adapted libraries (from Step 2.18)
- 80% Ethanol (from Step 2)
- Equilibrated DNA Purification Beads (from Step 2)
- Molecular biology grade water
- Optional 10 mM Tris-HCl pH 8 or Buffer EB
- Equinox Library Amp Mix (2x)
- From the Twist UMI Adapter System:
 - Twist UDI Primers

⚠ IMPORTANT: Use of Amplification Primers, ILMN tubes 100220, 100583 contained in the Twist Library Preparation EF Kit 1 are not required. Using these primers with the Twist UMI Adapter System will result in a failed PCR amplification.

Before You Begin

Thaw on ice:

- Twist UDI Primers (plate with single use primers)
- Equinox Library Amp Mix (2x)

PREPARE THE THERMAL CYCLER AND PCR MIX

3.1

Program a thermal cycler with the following conditions. Set the temperature of the heated lid to 105°C.

STEP	TEMPERATURE	TIME	NUMBER CYCLES
1 Initialization	98°C	45 seconds	1
2 Denaturation Annealing Extension	98°C	15 seconds	6–8*
	60°C	30 seconds	
	72°C	30 seconds	
3 Final Extension	72°C	1 minute	1
4 Final Hold	4°C	HOLD	—

*6–8 cycles is recommended when starting with 50–100 ng of high quality gDNA.

**PERFORM PCR**

3.2 Add 10 µl of Twist UDI Primer from the provided 96-well plate to each of the gDNA libraries from Step 2.18 and mix well by gentle pipetting.

NOTE: For index selection and multiplexing refer to the pooling guidelines in the Appendix.

3.3 Add 25 µl of Equinox Library Amp Mix (2x) to the gDNA libraries from Step 3.2 and mix well by gentle pipetting.

3.4 Pulse-spin sample plate or tube and immediately transfer to the thermal cycler. Start the program.

3.5 Remove the sample(s) from the block when the thermal cycler program is complete. Proceed to purification.

PURIFY

3.6 Vortex the pre-equilibrated DNA Purification Beads until mixed.

3.7 Add 50 µl (1x) of homogenized DNA Purification Beads to each ligation sample from Step 3.5. Mix well by vortexing.

3.8 Incubate the samples for 5 minutes at room temperature.

3.9 Place the samples on a magnetic plate for 1 minute.

3.10 The DNA Purification Beads form a pellet, leaving a clear supernatant. Without removing plate or tubes from the magnetic plate, remove and discard the supernatant.

3.11 Wash the bead pellet by gently adding 200 µl freshly prepared 80% ethanol (do not disturb the pellet), incubate for 1 minute, then remove and discard the ethanol.

3.12 Repeat this wash once, for a total of two washes, while keeping the samples on the magnetic plate.

3.13 Carefully remove all remaining ethanol with a 10-µl pipet, making sure not to disturb the bead pellet.

NOTE: Before pipetting, the bead pellet may be briefly spun to collect ethanol at the bottom of the plate or tube and returned to the magnetic plate.

3.14 Air-dry the bead pellet on the magnetic plate for 5 minutes or until the bead pellet is dry. Do not overdry the bead pellet.

3.15 Remove the plate or tubes from the magnetic plate and add 22 µl water, 10 mM Tris-HCl pH 8, or Buffer EB to each sample. Mix by pipetting until homogenized.


3.16 Incubate at room temperature for 2 minutes.

3.17 Place the plate or tubes on a magnetic plate and let stand for 3 minutes or until the beads form a pellet.

3.18 Transfer 20 μl of the clear supernatant containing the Amplified Indexed Libraries to a clean thin-walled PCR 0.2-ml strip-tube or well of a 96-well thermal cycling plate, making sure not to disturb the bead pellet.

PERFORM QC

3.19 Quantify and validate the size range of each library using the Thermo Fisher Scientific Qubit dsDNA Broad Range Quantitation Assay and Agilent DNA 7500 Assay. Final concentration values should be $\geq 80 \text{ ng}/\mu\text{l}$, and average fragment length should be 375–425 bp using a range setting of 150–1000 bp.

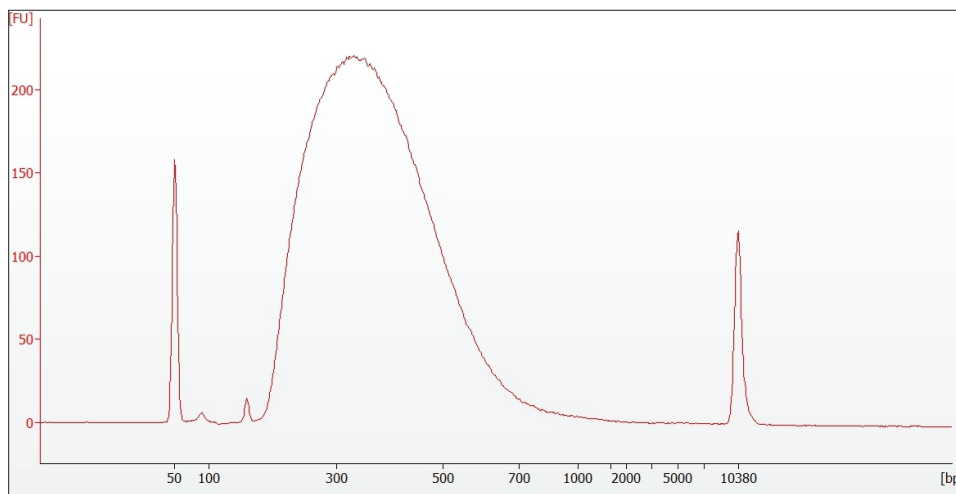

Electropherogram generated by an Agilent 7500 DNA analysis of gDNA library samples that were prepared as described. Note the single prominent peak at ~400 bp.

NOTES: If the average fragment length is not in the range of 375–425 bp:

- Input DNA concentration may be inaccurate. Typically, using too much DNA leads to shorter fragments, while not using enough DNA leads to larger fragments.
- The presence of cations and chelators may also affect the average fragment length.
- If neither of the above factors apply, optimize the 32°C fragmentation in Step 1.1 by changing the time in 3 minute increments—increase time to produce shorter fragments, and decrease time to produce longer fragments.

You can proceed with concentrations lower than 80 $\text{ng}/\mu\text{l}$, but low concentrations may reflect inefficient sample preparation and can result in low library diversity after hybridization.

STOPPING POINT: If not proceeding immediately to a Twist Target Enrichment System, store the amplified indexed libraries at -20°C .

END OF WORKFLOW



APPENDIX

UDI SEQUENCES

For a complete guide of the Twist UDI sequences please refer to PDF document DOC-001129 or Excel file DOC-001130. Both files are available for download at twistbioscience.com/resources.

POOLING GUIDELINES

Twist UDI primers are base balanced for next generation sequencing on a column basis. When pooling unique dual-indexed libraries for 8-plex hybridization, it is recommended that libraries be selected from a single column. Multiple columns may be selected in any desired combination across a single plate or multiple plates for sequencing.

Table 1. Twist UDI primer plate layouts and pooling guidelines.

Twist UMI Adapter System: TruSeq Compatible, 16 Samples (105040)

	1	2	3	4	5	6	7	8	9	10	11	12
A	1	9										
B	2	10										
C	3	11										
D	4	12										
E	5	13										
F	6	14										
G	7	15										
H	8	16										

*PLEASE NOTE: The indexes in the 16 sample plate are not the same in 96 samples, Plate A.



APPENDIX

Twist UMI Adapter System: TruSeq Compatible, 96 Samples, Plates A to D (105041, 105042, 105043, 105044)

Plate A.

	1	2	3	4	5	6	7	8	9	10	11	12
A	1	9	17	25	33	41	49	57	65	73	81	89
B	2	10	18	26	34	42	50	58	66	74	82	90
C	3	11	19	27	35	43	51	59	67	75	83	91
D	4	12	20	28	36	44	52	60	68	76	84	92
E	5	13	21	29	37	45	53	61	69	77	85	93
F	6	14	22	30	38	46	54	62	70	78	86	94
G	7	15	23	31	39	47	55	63	71	79	87	95
H	8	16	24	32	40	48	56	64	72	80	88	96

Plate B.

	1	2	3	4	5	6	7	8	9	10	11	12
A	97	105	113	121	129	137	145	153	161	169	177	185
B	98	106	114	122	130	138	146	154	162	170	178	186
C	99	107	115	123	131	139	147	155	163	171	179	187
D	100	108	116	124	132	140	148	156	164	172	180	188
E	101	109	117	125	133	141	149	157	165	173	181	189
F	102	110	118	126	134	142	150	158	166	174	182	190
G	103	111	119	127	135	143	151	159	167	175	183	191
H	104	112	120	128	136	144	152	160	168	176	184	192

Plate C.

	1	2	3	4	5	6	7	8	9	10	11	12
A	193	201	209	217	225	233	241	249	257	265	273	281
B	194	202	210	218	226	234	242	250	258	266	274	282
C	195	203	211	219	227	235	243	251	259	267	275	283
D	196	204	212	220	228	236	244	252	260	268	276	284
E	197	205	213	221	229	237	245	253	261	269	277	285
F	198	206	214	222	230	238	246	254	262	270	278	286
G	199	207	215	223	231	239	247	255	263	271	279	287
H	200	208	216	224	232	240	248	256	264	272	280	288

Plate D.

	1	2	3	4	5	6	7	8	9	10	11	12
A	289	297	305	313	321	329	337	345	353	361	369	377
B	290	298	306	314	322	330	338	346	354	362	370	378
C	291	299	307	315	323	331	339	347	355	363	371	379
D	292	300	308	316	324	332	340	348	356	364	372	380
E	293	301	309	317	325	333	341	349	357	365	373	381
F	294	302	310	318	326	334	342	350	358	366	374	382
G	295	303	311	319	327	335	343	351	359	367	375	383
H	296	304	312	320	328	336	344	352	360	368	376	384

END OF APPENDIX